



Walk About™ Insulative Scoop Tray

Cat. No. 39438-001 Walk About Insulative Scoop Tray with Beads
Cat. No. 39438-002 Walk About Insulative Scoop Tray

Description

The Walk About Scoop Tray is a revolutionary way to keep your samples at temperature as you work under the hood, at the bench, or as you travel between labs. It's a quarter the size of a traditional lab bucket which makes it easy to use in tight benchtop spaces. With its superior insulative properties, it keeps your samples and reagents at temperature after removing them from the bead bath, Chill Bucket, or the refrigerator.

Walk About™ Features

- ✓ One half to one liter capacity
- ✓ Expanded PVC construction
- ✓ Compatible from -196°C to 93°C
- ✓ Durable and lightweight
- ✓ Small size for use under the hood or on the bench
- ✓ Molded one-piece construction
- ✓ Accepts and supports any size and shape vessel including non-water-tight vessels
- ✓ Washes and disinfects easily
- ✓ Condensation free

Setup

Use the following procedures to set up your Walk About Scoop Tray.

For 37°C:

- Store empty Walk About at ambient temperature.
- Scoop approximately 0.5L of Lab Armor Beads from bead bath set to 37°C.
- Add samples.
- After use, disassemble the Walk About and return the Beads to the bead bath.

Alternative Method for 37°C:

- Store Walk About containing beads at 37°C in an oven or other heating device.
- Remove Walk About from heating device and add samples.
- After use, return Walk About to heating device.

For 4°C to 8°C use:


- Store Walk About containing Beads at 0°C-4°C in refrigeration unit.
- Remove Walk About from refrigeration unit and add samples.
- After use, return Walk About to refrigeration unit.

Tips and Hints

- ✓ Store Beads at the desired working temperature.
- ✓ If necessary, wash Beads and tray with soap and water. **Completely** dry Beads before returning to the tray.
- ✓ Avoid using strong acids, bases, including bleach solutions, and detergents.

If you have any doubts about which cleaning products to use, we recommend our Lab Armor® Cleaning Products.

Part ID	Description	Part ID	Description
1200518	Lab Armor® Bead Soap Wash	1200520	Lab Armor® Cleanser

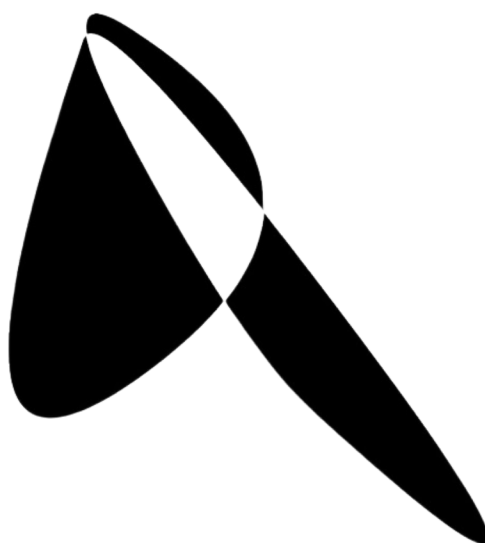

Technical Support

For additional product and technical information, such as product manuals or technical articles on the use of the Walk About Tray and Lab Armor Beads for common laboratory applications, please visit our website at www.labarmor.com. For further assistance, please email us at labarmor@sheldonmfg.com.

www.labarmor.com | labarmor@sheldonmfg.com | +1 (503) 640-3000 | +1 (800) 322-4897 (U.S. Toll-Free)

Notifications

The trademarks mentioned herein are the property of Lab Armor, LLC. Lab Armor® Beads are patented. For laboratory use only. Not for personal use. © 2010-2025 Lab Armor. All rights reserved.