



## Stay Temp™ Tray

Cat. No. 39956-050     1L Stay Temp Tray  
 Cat. No. 39956-001     2L Stay Temp Tray  
 Cat. No. 39956-003     4L Stay Temp Tray

### Description

Stay Temp trays are an upgrade that keeps samples and reagents in your refrigerator, incubator, and laboratory ovens at temperature despite normal hourly fluctuations in air temperature. Samples will be protected from defrost cycles and constant door opening and closing. StayTemp trays also eliminate the need for racks.

### Stay Temp™ Features

- ✓ Stainless steel construction
- ✓ Keeps samples organized
- ✓ For use in refrigerators, incubators, & ovens
- ✓ Maintains cold or heat with high efficiency
- ✓ Accepts and supports any size and shape vessel, including non-water-tight vessels
- ✓ Hand wash, autoclave, towel dry
- ✓ 3 sizes available, does NOT include lid

### Setup

Use the following procedures to set up your Stay Temp system.

1. Place StayTemp tray on workbench or counter.
2. Cut the corner of one side of the Lab Armor Bead Bag and pour beads into tray.
3. Place filled tray into desired equipment and allow to equilibrate 4-6 hours.
4. Insert vessels into Beads until firmly supported.

### Tips and Hints

- ✓ Store Beads at the desired working temperature.
- ✓ If necessary, wash Beads and tray with soap and water. **Completely** dry Beads before returning to the tray.
- ✓ Avoid using strong acids, bases, including bleach solutions, and detergents.

*If you have any doubts about which cleaning products to use, we recommend our Lab Armor® Cleaning Products.*

Part ID	Description	Part ID	Description
1200518	Lab Armor® Bead Soap Wash	1200520	Lab Armor® Cleanser

### Technical Support

For additional product and technical information, such as product manuals or technical articles on the use of the Stay Temp Tray and Lab Armor Beads for common laboratory applications, please visit our website at [www.labarmor.com](http://www.labarmor.com). For further assistance, please email us at [labarmor@sheldonmfg.com](mailto:labarmor@sheldonmfg.com).

[www.labarmor.com](http://www.labarmor.com) | [labarmor@sheldonmfg.com](mailto:labarmor@sheldonmfg.com) | +1 (503) 640-3000 | +1 (800) 322-4897 (U.S. Toll-Free)

### Notifications

The trademarks mentioned herein are the property of Lab Armor, LLC. Lab Armor® Beads are patented. For laboratory use only. Not for personal use. © 2010-2025 Lab Armor. All rights reserved.