



Lot Number:

StayTemp™ Equipment Tray

Cat. No. 39956-050

Cat. No. 39956-001

Cat. No. 39956-003

1L StayTemp™ Equipment Tray

2L StayTemp™ Equipment Tray

4L StayTemp™ Equipment Tray

Description

StayTemp™ Equipment trays are an upgrade that keeps samples and reagents in your refrigerator, incubator, and laboratory ovens at temperature despite normal hourly fluctuations in air temperature. Samples will be protected from defrost cycles and constant door opening and closing. StayTemp™ equipment trays also eliminate the need for racks.

StayTemp™ Equipment Tray features:

- ✓ Stainless Steel construction
- ✓ Keeps samples organized
- ✓ For use in refrigerators, incubators, & ovens
- ✓ Maintains cold or heat with high efficiency
- ✓ Accepts and supports any size and shape vessel including non-water-tight vessels
- ✓ Hand wash, autoclave, towel dry
- ✓ 3 sizes available, does not include lid

Tips and Hints

- ✓ If necessary, wash Beads and tray with soap and water. *Completely dry* Beads before returning to the tray
- ✓ Avoid using strong acids, bases, including bleach solutions, and detergents
- ✓ For more info, visit <http://www.support.labarmor.com/products.html>

Setup

Use the following procedures to set up your StayTemp™ system.

- ✓ Place StayTemp™ tray on workbench or counter.
- ✓ Cut the corner off one side of the Lab Armor® Bead Bag and pour beads into tray
- ✓ Place filled tray into desired equipment and allow to equilibrate 4-6 hours
- ✓ Insert vessels into Beads until firmly supported

Notifications

Lab Armor® Beads are Patent Pending. For laboratory use only. Not for personal use. © 2011 Lab Armor®. All Rights Reserved. Version 2.3 REV: 042810